

EFS pacing kit for FDSS/μCELL

Specialized for Cardiomyocyte Assays


Overview

- Cardiomyocyte Pacing
- Electric Stimulation of Human iPSC-derived Neurons

Electric Field Stimulation (EFS) pacing kit for FDSS/μCELL is an option to have an ability to stimulate cardiomyocytes electrically by the 96-electrode array to pace beatings of the cells, in which the cells are cultivated in a 96-well plate format. This option can be associated with other option available in FDSS/μCELL, such as assay plate heating unit and fast data acquisition software (up to 120 Hz frame reading), each of which could make more accurate monitoring of intracellular molecular events such as Ca²⁺ transients in cardiomyocytes. The EFS option with FDSS/μCELL should be useful in the drug discovery and in cardiomyocyte research. The EFS system also can stimulate neurons electrically.

Application 1

- Cardiomyocyte Pacing
Measurements of intracellular Ca²⁺ concentration changes

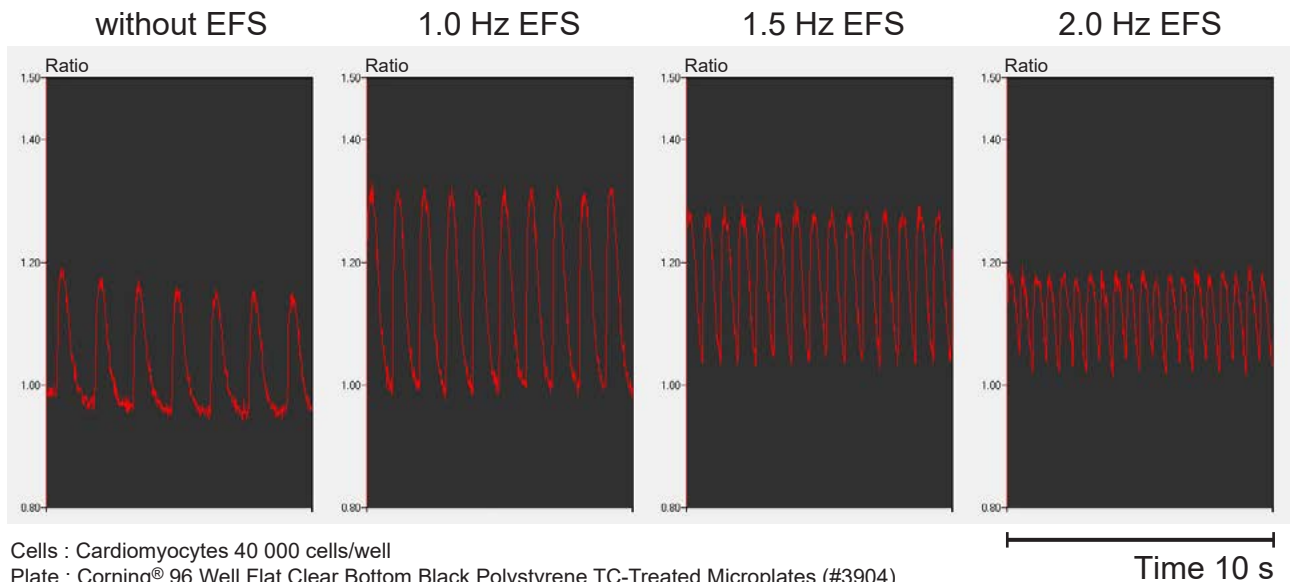

Cells : Cardiomyocytes 40 000 cells/well

Plate : Corning[®] 96 Well Flat Clear Bottom Black Polystyrene TC-Treated Microplates (#3904)

Dye : Cal-520/AM final 2 μM, Probenecid final 1.25 mM, Loading 45 min

System : FDSS/μCELL, EM-CCD 2×2 binning, exposure time 16 ms

EFS (Electric Field Stimulation) : 5 V, 10 ms Interval, 1.0/1.5/2.0 Hz, 60/90/120 times

Caution Notice:

The FDSS/μCELL EFS system should not be used for optically detecting/monitoring change in transmembrane potential of the cells.

The FDSS/μCELL EFS system should not be used on any cell or cells in which the user or anyone else has expressed target ion channels.

FDSS Option

Application 2

● Electric Stimulation of Human iPSC-derived Neurons Measurements of intracellular Ca²⁺ concentration changes


Cells : Peri.4U® (human iPSC-derived Peripheral neurons, Ncardia) 40 000 cells/well

Plate : Corning 96 Well Flat Clear Bottom Black Polystyrene TC-Treated Microplates (#3904)

Dye : Cal-520/AM final 2 μM, Probenecid final 1.25 mM, Loading 45 min

System : FDSS/μCELL, EM-CCD 1×1 binning, exposure time 153 ms

EFS (Electric Field Stimulation) : 20 V, 1 ms Interval, 10/20/30/40 Hz, 50 times at each stimulation

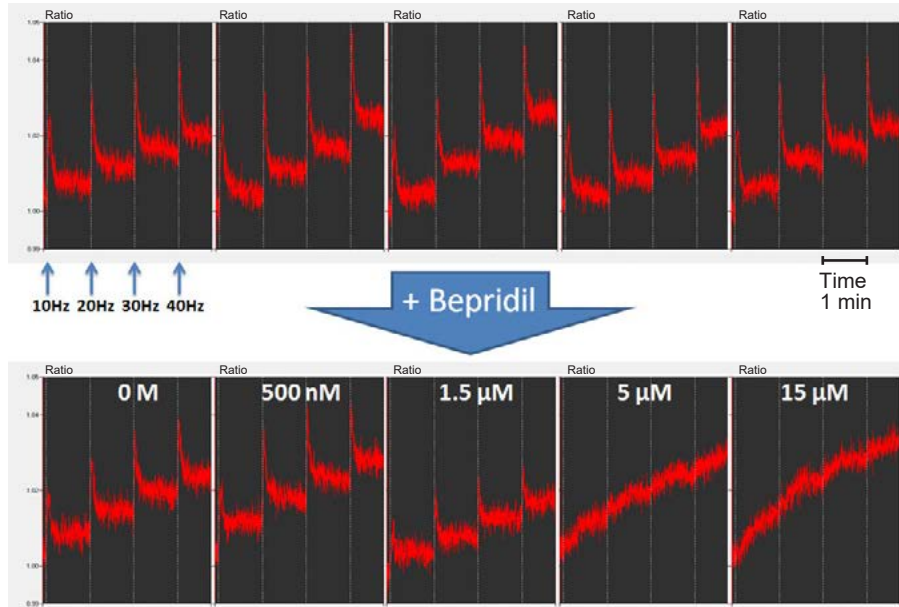

Electric stimulations were added to human iPSC-derived peripheral neurons at 10, 20, 30, and 40 Hz sequentially. At concentration more than 5 μM Bepridil, the intracellular Ca²⁺ concentration changes were completely inhibited.

▲ FDSS/μCELL

- Fluorescence / Luminescence measurements while electric field stimulation.
- Variable EFS frequency.
- Variable EFS voltage.

* EFS pacing kit option requires fluorescence / luminescence sensor and the latest data analysis computer.
Please contact your local sales representative for more detail.

| Part number | Description |
|-------------|---|
| A11529-14 | Heater unit |
| U8524-11 | FDSS Software option High Speed Acquisition Option |
| U8524-12 | FDSS Software option Waveform Analysis software for cardiomyocyte |
| M13040-01 | EFS pacing system |

★ FDSS is a registered trademark of Hamamatsu Photonics K.K. (EU, Japan, U.S.A)

★ Hamamatsu Photonics K.K. and its subsidiaries or representatives do not warrant that the use of our products and those goods does not infringe the intellectual property rights of third parties.

★ Product and software package names noted in this documentation are trademarks or registered trademarks of their respective manufacturers.

★ Subject to local technical requirements and regulations. Availability of products included in this promotional material may vary. Please consult with your local sales representative.

● Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions.

● Specifications and external appearance are subject to change without notice.

© 2020 Hamamatsu Photonics K.K.

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

Systems Division

812 Joko-cho, Higashi-ku, Hamamatsu City, 431-3196, Japan, Telephone: (81)53-431-0124, Fax: (81)53-433-8031, E-mail: export@sys.hpk.co.jp

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, Bridgewater, NJ 08807, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-265-8 E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: infos@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire AL7 1BW, UK, Telephone: (44)1707-294888, Fax: (44)1707-325777 E-mail: info@hamamatsu.co.uk

North Europe: Hamamatsu Photonics Norden AB: Torshamnsgatan 35 16440 Kista, Sweden, Telephone: (46)8-509 031 00, Fax: (46)8-509 031 01 E-mail: info@hamamatsu.se

Italy: Hamamatsu Photonics Italia S.r.l.: Strada della Moia, 1 int. 6, 20020 Arese (Milano), Italy, Telephone: (39)02-93 58 17 33, Fax: (39)02-93 58 17 41 E-mail: info@hamamatsu.it

China: Hamamatsu Photonics (China) Co., Ltd.: 1201 Tower B, Jiaming Center, 27 Dongsanhuan Beilu, Chaoyang District, 100020 Beijing, P.R.China, Telephone: (86)10-6586-6006, Fax: (86)10-6586-2866 E-mail: hpc@hamamatsu.com.cn

Taiwan: Hamamatsu Photonics Taiwan Co., Ltd.: 8F-3, No.158, Section2, Gongdao 5th Road, East District, Hsinchu, 300, Taiwan R.O.C. Telephone: (886)3-659-0080, Fax: (886)3-659-0081 E-mail: info@hamamatsu.com.tw

Cat. No. SBIS0114E03
MAR/2020 HPK
Created in Japan